

Free SO₂

Cat. no. 301

Intended Use

AV-Free SO₂ is intended for measuring the level of unbound Sulfur Dioxide in wine, grape juice and must.

Methodology

AV-Free SO₂ is based on the reduction in color exhibited by a dye when it reacts with sulfur dioxide.



Sample

Samples of wine and must should be used as is. The ACCUVIN AV-Free SO₂ test minimizes the usual interferences from colored and turbid samples. Samples do not have to be pre-filtered or treated with color removing substances such as activated carbon or polyamide powder. Sample temperature may be from 0°C - 35°C (32°F - 95°F).

Procedure

Note: The ACCUVIN AV-Free SO₂ test covers two ranges. A high range test (tubes with red caps) covers the range of 40 – 130 ppm, and is usually applicable for sweet wines and for wines with a pH of 3.7 or higher. The low range test (tubes with green caps) covers the range 0 – 40 ppm, and is usually applicable for wines with a pH of 3.6 or less.

1. Pour clear liquid from tube with black cap into powder into a tube with a green cap or a tube with a red cap.
2. Replace green or red cap and shake tube until all powder is dissolved, about 1 minute.
3. Squeeze upper sampler bulb. Dip sampler tip into grape juice, must or wine sample, then **slowly release bulb** to aspirate sample. **Wipe sampler tip to remove excess droplets.** (If you prefer to use an air displacement pipette, set sample volume at 91 µL.)
4. Open sample tube. Transfer sample to test tube by placing sampler tip into the test reagent and squeezing sample bulb only once. Withdraw sampler prior to releasing sampler bulb. Note that only sample present in the sampler tip will be dispensed. Replace sample tube cap. Shake tube. Wait 5 min. for color development.
5. Determine sample Free SO₂ level in ppm by comparing the developed color to the color chart on the test strip container. Read tube color by holding tube approximately 1 inch (2.5 cm.) above a white background. If test strip color falls between two color chips select an intermediate value for the sample Free SO₂ level. (Since fluorescent lights have a green cast, color matching is best under incandescent or natural lighting.)

Storage

Store away from direct sunlight at temperatures below 80°F. Product is satisfactory until the date printed on the test tube container label.

www.ACCUVIN.com

for technical inquiries: email: contact@accuvin.com

Seller's limitation of liabilities: Every effort has been made to ensure the material contained in this informational insert and the results obtained with AV test strips are as accurate as possible, but no warranty or fitness is implied. Buyer shall not in any event be entitled to, and seller shall not be liable for, indirect, special, incidental or consequential damages of any nature including, without being limited to, loss of profit, promotional or manufacturing expenses, overhead, injury to reputation or loss of customers. Buyer's recovery from seller for any claim shall not exceed buyer's purchase price for the products irrespective of the nature of the claim, whether in contract, tort, warranty, or otherwise.

Interpretation

Sulfur dioxide is a chemical not normally present in grapes, but when present it can have a powerful antibacterial effect.

One source of sulfur dioxide in musts is from yeasts during alcoholic fermentation. Although not a prevalent problem, the levels produced in this manner may be sufficient to cause inhibition of malolactic bacteria. These levels are dependent on the yeast strain, the availability of nutrients, and the presence of compounds in the must to which sulfur dioxide binds.^{1, 2, 5}

A more common source of sulfur dioxide in musts and wines is through addition by the winemaker. These additions are made for the following reasons:

- At crush/destemming to inhibit native bacteria, especially lactic acid bacteria, which could cause formation of undesirable compounds such as acetic acid.
- At crush/destemming to inhibit browning, especially in white grapes
- After alcoholic fermentation to inhibit lactic acid bacteria and reduce the likelihood of unplanned malolactic fermentation by malolactic organisms.
- During aging and just prior to bottling to reduce enzymatic and chemical oxidation and control bacteria.

The total amount of sulfur dioxide present in must and wine is not all useful as a preservative. Some is bound to other chemical components, reducing its effectiveness as a preservative by 30 – 70%. Also, the degree of binding can change over time, affecting the ability of an adequate dose to continue providing its preservative properties. The effectiveness of sulfur dioxide as a preservative is also pH-dependent, with low pH wines requiring lower levels, and high pH wines requiring significantly greater concentrations to achieve the same effectiveness as in a typical wine with a pH of 3.3 - 3.4.

Summary Interpretation for Most Wines

(Because of varietal & stylistic differences, growers & winemakers should make final interpretations.)

Free SO₂	Interpretation for wines at pH 3.3 – 3.5
20 - 30 ppm	Level to control harmful microorganisms in fresh musts for red wine from clean, sound fruit pre-fermentation w/o inhibiting MLF ⁴
30 - 50 ppm	Level to control browning in white musts, harmful microorganisms in fresh juice from clean, sound fruit ³
70 - 100 ppm	Level to control browning in white musts, harmful microorganisms in fresh juice from fruit showing evidence of spoilage or rot ^{3,4}
20 – 40 ppm	Level desired for bottling ²

Application Table

pH	% molecular SO ₂	free SO ₂ needed to achieve molecular SO ₂ of:		
		<u>0.5 ppm</u>	<u>0.8 ppm</u>	<u>2.0 ppm*</u>
2.9	7.5	7	11	27
3.0	6.1	8	13	33
3.1	4.9	10	16	41
3.2	3.9	13	21	51
3.3	3.1	16	26	64
3.4	2.5	20	32	80
3.5	2.0	25	40	100
3.6	1.6	31	50	125
3.7	1.3	39	63	157
3.8	1.0	49	79	197
3.9	0.8	62	99	248
4.0	0.7	78	125	312

* for sweet white wines

References

1. Margalit, Y., Concepts in Wine Chemistry, The Wine Appreciation Guild, San Francisco, **1997**. p. 18
2. Zoecklein, B.W., Fugelsang, K.C., Gump, B.H. and Nury, F.S., *Wine Analysis and Production*, Chapman and Hall, New York, **1995**. pp. 192 - 196.
3. Dharmadhikari, M., "Vineyard and Vintage View," 14(3) 2 **1999**
4. Dharmadhikari, M., "Vineyard and Vintage View," 15(3) 1 **2000**
5. Henick-Kling, T., and Y. H. Park, "Considerations for the use of yeast and bacterial starter cultures: SO₂ and timing of inoculation," Am. J. Enol. Vitic., 45 (4): 464 – 469 **1994**.